

 Phase: **MEGA-DEX DET Beta**
Catalog Code: DEX-DET-B-025-025-25

 Phase: **MEGA-DEX DMT Beta**
Catalog Code: DEX-DMT-B-025-025-25

 Phase: **MEGA-DEX DAC Beta**
Catalog Code: DEX-DAC-B-025-025-25

Columns Dimensions: **0.25mm, 0.25µm, 25m.**

Camphor


Chiral Separation on MEGA-DEX Columns

Conditions:

Injection: Split 250°C, 0.5µL, 1:50 Split Ratio.

Sample dilution: 1mg/1mL in n-Hexane.

Contact us for additional information.

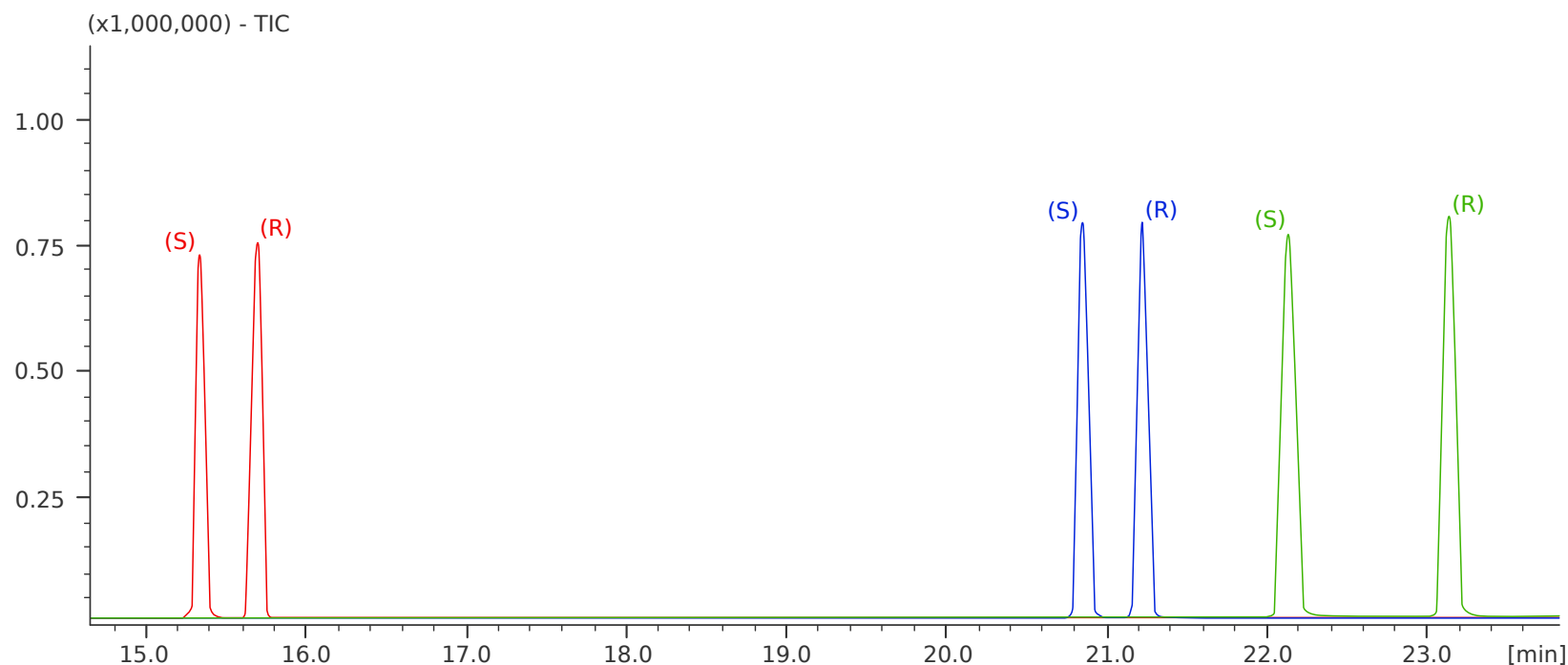

Acknowledgement: Prof. C.Bicchi, C.Cordero et al. - Università di Torino, Dipartimento di Scienza e Tecnologia del Farmaco, Via Pietro Giuria, 9 - Torino - Italy.

