

Residual Solvents from Food Packaging

Column: **MEGA-VOC 2 FAST - 0.20mm, 1.00µm, 10m**

Catalog Code: F-VOC2-020-100-10

Conditions:

Injection: Split 250°C.

Detector: FID 250°C.

Oven Temp.: 35°C (0.5min), 15°C/min, 100°C.

Carrier Gas: Hydrogen, 50 kPa.

Identificazione:

1. Methanol
2. Ethanol
3. Isopropanol
4. Methyl Ethyl Ketone (MEK)
5. Ethyl Acetate
6. Isobutanol
7. Cyclohexane
8. 1-Methoxy-2-Propanol
9. Ethoxy Propanol
10. Isobutyl Acetate
11. N-Butyl Acetate

Analysis carried out with DANI MASTER-GC.

